Clean-Trace™ Test


-  **(EN)** Surface Positive Control
-  **(FR)** Témoin positif pour les surfaces
-  **(DE)** Oberflächen-Positivkontrolle
-  **(IT)** Controllo positivo delle superfici
-  **(ES)** Control positivo para superficies
-  **(NL)** Oppervlakpositieve controle
-  **(SV)** Positivt yt-test
-  **(DA)** Overflade Positiv Kontrol
-  **(NO)** Positiv kontroll til overflate
-  **(FI)** Surface positiivinen kontrolli
-  **(PT)** Controle Positivo de Superfície
-  **(EL)** Θετικής Επιβεβαίωσης για Επιφάνειες
-  **(PL)** Kontrola dodatnia powierzchni
-  **(HU)** Felszíni pozitív kontroll
-  **(CS)** Povrchová pozitivní kontrola
-  **(RO)** Control pozitiv al suprafețelor
-  **(RU)** Положительный контроль поверхности
-  **(TR)** Yüzey Pozitif Kontrolü
-  **(JA)** 表面ポジティブコントロール
-  **(ZH)** 表面采样棒阳性对照

Clean-Trace™ Test

Product Instructions

Surface Positive Control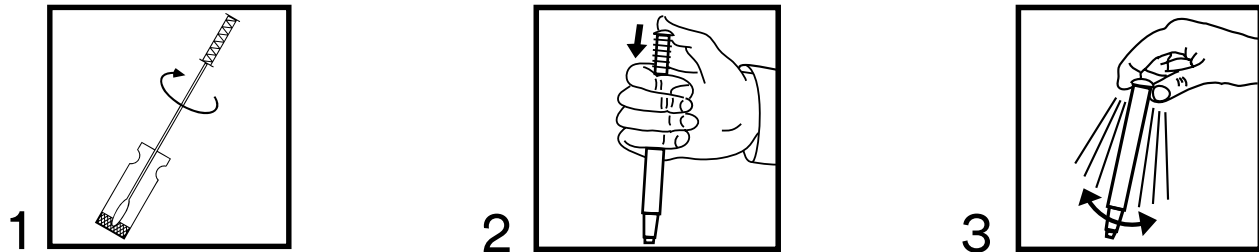

Product Description and Intended Use

The measurement of the standard concentration of Adenosine Triphosphate (ATP) or protein by 3M™ Clean-Trace™ Surface Tests is used as an assurance of reagent and in the case of, the ATP tests which require a 3M™ Clean-Trace™ Luminometer, instrument function. The test kit contains 10 ready to use reagent vials. Each vial contains 3.0 ng or 5×10^{-12} moles of ATP and 1 mg of protein (Bovine Serum Albumin). A pre-moistened Test Swab is used to swab the ATP/protein pellet and interior of the 3M™ Clean-Trace™ Surface Positive Control vial. 3M Clean-Trace Surface Test devices are then activated and measured as per the test device instructions. The signal obtained should be above a specified level, (see below) which indicates that both the reagents and (in the case of the ATP tests) instrument are operating correctly. The unique format of this product provides the user with an ideal and convenient method of assuring that all components used during testing are operating correctly, and that the operator is using the correct method. 3M xClean-Trace Surface Positive Control can be used in conjunction with the 3M™ Clean-Trace™ Surface ATP Test, 3M™ Clean-Trace™ Surface Protein Plus Test, 3M™ Clean-Trace™ Surface Protein High Sensitivity Test, and 3M™ Clean-Trace™ Surface Protein Allergen Test to rapidly assess that both reagents and instruments (if required) are operating satisfactorily and have been stored and used correctly.

⚠ Warnings

Consult the Safety Data Sheet for additional information and local regulations for disposal. ATP is a common substance. It is present on skin, hair and on many surfaces. Do not touch the swab or the interior surfaces of the positive control vials. Do not use devices past expiry date.

Notice

To reduce the risk of inaccurate results 3M Clean-Trace Surface Positive Control Kit should be stored at 2°-8°C (36°-46°F). To avoid inaccurate results, keep reconstituted 3M Clean-Trace Surface Positive Control Kit at refrigerated temperatures 2°-8°C (36°-46°F). To reduce the risks associated with cross contamination while preparing the assay the use of gloves is recommended. To reduce the risk of inaccurate results do not shake the reconstituted reagents. Follow instructions for reconstituting reagents.

User Responsibility

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.3M.com/foodsafety, or contact your local 3M representative or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any 3M Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

Limitation of Warranties / Limited Remedy

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, 3M DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any 3M Food Safety Product is defective, 3M or its authorized distributor will, at its option, replace, or refund the purchase price of the product. These are your exclusive remedies. You must promptly notify 3M within sixty days of discovery of any suspected defects in a product and return it to 3M. Please all Customer Service (1-800-328-1617 in the U.S.) or your official 3M Food Safety representative for a Returned Goods Authorization.

Limitation of 3M Liability

3M WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall 3M's liability under any legal theory exceed the purchase price of the product alleged to be defective.

Storage and Disposal

For maximum shelf life, store between 2°-8°C (36°-46°F). Expiry date and lot number are noted on each package. After use, may contain microorganisms that may be a potential biohazard. Follow current industry standards for disposal. Do not freeze. Protect from direct sunlight.

Instructions for Use

Before the 3M Clean-Trace Surface Positive Control is used, it is important to ensure your 3M Clean-Trace Luminometer is switched on and initialized. Refer to the manual provided with the instrument for full details.

1. Place the 3M Clean-Trace Test devices and positive control vials at room temperature for at least 10 minutes before use.
2. Gently tap the positive control vial to ensure that the reagent pellet is at the bottom.
3. Carefully remove the stopper from the vial, avoid contaminating the opening to the vial.
4. Immediately before use, remove the swab test device from the foil pouch. Grip the sample stick handle and remove the swab from the device.
5. For the 3M Clean-Trace Surface Protein Plus Test, 3M Clean-Trace Surface Protein High Sensitivity Test and 3M Clean-Trace Surface Protein Allergen Test, pre-moisten the swab head as per test device instructions.
6. Swab around the interior of the vial, twisting the swab to pick up the entire pellet [figure 1].
7. Reinsert the sample stick into the device with the handle inserted to the original position of the unused device.
8. To process the sample, push down firmly on the top of the sample stick handle [figure 2]. The handle will slide into the test device tube and the top of the handle should be level with the top of the test device tube when fully depressed. Grip the top of the device and shake rapidly side-to-side for at least five seconds to mix the sample and reagent. [figure 3]
9. Read the swab test device as per the swab product instructions.
10. Replace the positive control vial stopper and dispose of the vial.

Interpretation

Testing using the 3M Clean-Trace Surface ATP Test should give readings greater than 10,000 RLU.

Results of testing using the 3M Clean-Trace Surface Protein Tests should give a dark purple color change within 10 minutes of activation with the exception of 3M™ Clean-Trace™ Surface Protein Instant (CIT50), which is not compatible with the 3M Clean-Trace Surface Positive Control.

If the tests fail to reach these levels, see the troubleshooting guide below and repeat the test procedure using a fresh vial of positive control reagent.

If low results are again achieved it may indicate a possible problem with either the reagent or instrument (if applicable).

Trouble Shooting Guide

The following section is a list of potential problems that may cause a low signal:

- Incomplete mixing - shake the device and read again
- Reagents have passed their expiration date
- Reagents stored incorrectly
- Reagents not allowed to reach room temperature before testing
- Poor swabbing technique - check that there is no residual pellet left in the bottom of the vial
- Contaminated luminometer
- Damaged luminometer

For further information, contact 3M Food Safety Technical Services.

Explanation of Symbols

www.3M.com/foodsafety/symbols

3M Food Safety

3M United States

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-800-328-6553

3M Canada

Post Office Box 5757
London, Ontario N6A 4T1
Canada
1-800-563-2921

3M Latin America

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-954-340-8263

3M Europe and MEA

3M Deutschland GmbH
Carl-Schurz-Strasse 1
D41453 Neuss/Germany
+49-2131-14-3000

3M United Kingdom PLC

Morley Street, Loughborough
Leicestershire
LE11 1EP
United Kingdom
+(44) 1509 611 611

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Euro Plaza
Gebaude J, A-1120 Wien
Kranichberggasse 4
Austria
+(43) 1 86 686-0

3M Asia Pacific

No 1, Yishun Avenue 7
Singapore, 768923
65-64508869

3M Japan

3M Health Care Limited
6-7-29, Kita-Shinagawa
Shinagawa-ku, Tokyo
141-8684 Japan
81-570-011-321

3M Australia

Bldg A, 1 Rivett Road
North Ryde, NSW 2113
Australia
61 1300 363 878

3M Health Care

2510 Conway Ave
St. Paul, MN 55144 USA
www.3M.com/foodsafety

© 2018, 3M. All rights reserved.
3M and Clean-Trace are trademarks of 3M. Used under license in Canada.
34-8723-5572-1

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3M Österreich GmbH